

HSM

Manual Vibratory Disc Mill



HERZOG

Efficient and reproducible fine grinding

Manual Vibratory Disc Mill Model HSM


The HSM Vibratory Disc Mill is the optimal machine for fine grinding of hard, medium-hard and brittle materials. The HSM provides highly efficient particle size reduction for a wide range of materials, including mineral materials, slags, ferroalloys, and cement products. The simple and ergonomic operability, the low noise level and small footprint make it the optimal machine for laboratory operation. At the same time, the machine can be used in industrial environments without any problems due to its extremely robust and resistant design.

Maximum flexibility


The operator can easily configure all grinding parameters on the control panel of the HSM and compile individual sample preparation programs. The programmable logic controller allows up to 30 different program sequences to be stored. The speed of the mill is continuously adjustable in a range between 750 and 1500 rpm. The automatic program sequence guarantees reproducible sample preparation and thus a high degree of accuracy and precision in material analysis. Due to the high rotation speed, even hard materials are ground down to the required particle size within short preparation times.

To enable analysis-neutral comminution of the sample material for different applications, grinding vessels and sets are offered in different material qualities. Optional versions are available in chrome steel, tungsten carbide, Colmonoy and agate.

Various grinding vessel sizes are available for efficient grinding of different sample quantities. In the chromium steel and tungsten carbide version, grinding vessel volumes of 10, 50, 100 and 250 ccm are available. In the Colmonoy version, grinding vessels with a volume of 100 or 250 ccm can be used. The appropriate grinding sets are adapted in such a way that an optimum grinding result is achieved with each grinding vessel size.

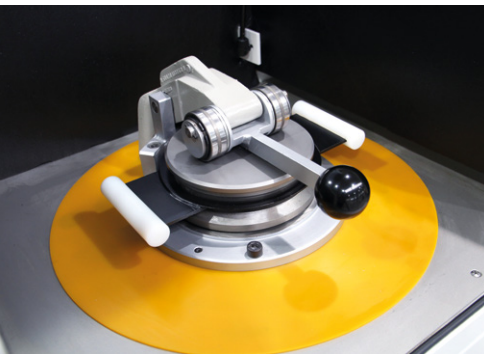


Manual clamping device

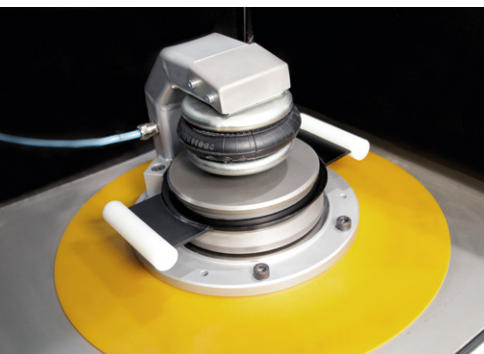

Pneumatic clamping device



Grinding vessel with carrying grips

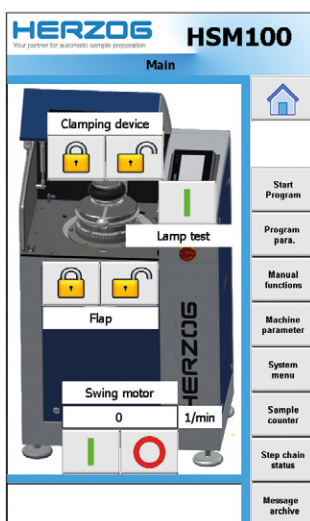

Easy operability due to clear menu navigation



Perfectly suited for a wide range of different applications

Great ergonomics and operability

A major focus was placed on the safe and ergonomic operation of the HSM. The working height has been adjusted to allow effortless loading and unloading of the grinding vessel. A broad carrying grip on both sides makes handling of the grinding vessel easier, resulting in greater comfort and safety during transport.

Accessibility to the machine has been further improved by enlarging the grinding chamber. The operator can open and close the machine door safely and comfortably with the help of gas springs. During machine operation, access to the machine chamber is prevented by a safety switch with guard locking. The completely enclosed grinding chamber and the sound-insulating lining of the housing result in a significant noise reduction.

In the basic HSM-H variant, the grinding vessel is clamped by means of an easy-to-operate quick-action clamping device. In the optional HSM-P variant, the grinding jar is automatically fixed by a pneumatic clamping device.

The machine is operated via the integrated touch-sensitive 7-inch color display. The panel size and clear menu navigation allow easy and intuitive access to all machine functions. The three buttons located on the front of the machine are programmable and allow the direct start of frequently used preparation programs.

Improved service properties

The HSM features improved maintenance and service characteristics. Optionally available machine feet enable the machines to be set up on uneven floors and allow easy adjustment to the operator's optimum working height.


The front- and side-mounted inspection hatches facilitate access to the interior of the machine, enabling operators to perform regular cleaning and inspection tasks with ease. In addition, any necessary service and repair work can be carried out quickly and efficiently.



Ergonomic and safe operation



Integrated 7" color panel

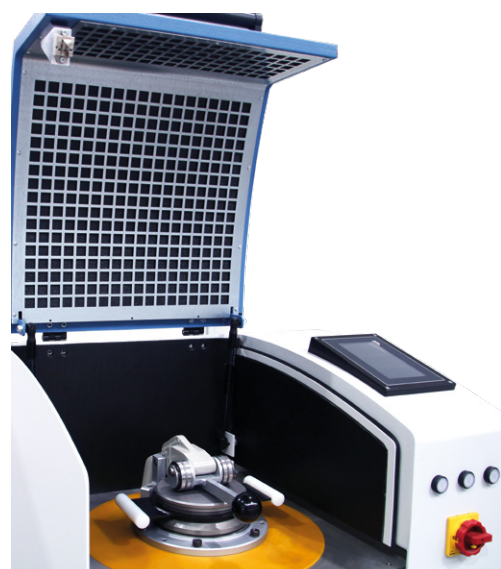

Optimal access to the grinding chamber

Technical data

Manual Vibratory Disc Mill HSM

Dimensions L x W x H

Machine	600 x 780 x 1275 mm
Machine including packing	1200 x 1200 x 1800 mm

Weight

Machine	275 kg
Machine including packing	413 kg

Power supply and consumption

Voltage	380–420 V, 50 Hz, 3-phase 380–480 V, 60 Hz, 3-phase 220–240 V, 50/60 Hz, 3-phase (Other voltages possible on request)
Center conductor	Not required
Power consumption	2.0 kVA

Electrical control cabinet (integrated)

PLC	SIMATIC S7-1200
Control voltage	24 V DC
Protection class	IP 54
Insulation class	B

Pneumatic supply and consumption (HSM-P only)

Pressure	Min. 5 bar, max. 10 bar
Consumption per sample	Approx. 10 dm ³

Machining parameters

Grinding time	1–900 S
Speed	750–1500 min ⁻¹ (infinitely variable)

Sample types

Material	Various hard, medium-hard and brittle materials such as minerals, slags, sinter, ferroalloys, cement, cement precursor and others, organic materials
Material hardness	Max. 9 Mohs
Input grain size	10 mm

Options

- Pneumatic clamping device (HSM-P)
- Grinding vessels and grinding set made of chrome steel (10, 50, 100 or 250 ccm)
- Grinding vessels and grinding set made of tungsten carbide (10, 50, 100 or 250 ccm)
- Grinding vessels and grinding set made of Colmonoy (100 or 250 ccm)
- Grinding vessel and grinding set made of agate (100 ccm)



The design of the machine complies with the applicable accident prevention and VDE regulations. We reserve the right to make technical modifications.

HSM/1.2021-D-1

HERZOG Maschinenfabrik GmbH & Co. KG

Auf dem Gehren 1
49086 Osnabrück
Germany
Phone +49 541 9 33 20
Fax +49 541 9 33 232

info@herzog-maschinenfabrik.de
www.herzog-maschinenfabrik.de

HERZOG Automation Corp.

16600 Sprague Road, Suite 400
Cleveland, Ohio 44130
USA
Phone +1 440 891 9777
Fax +1 440 891 9778

info@herzogautomation.com
www.herzogautomation.com

HERZOG Japan Co., Ltd.

3-7, Komagome 2-chome
Toshima-ku, Japan
Tokyo 170-0003
Phone +81 3 5907 1771
Fax +81 3 5907 1770

info@herzog.co.jp
www.herzog.co.jp

HERZOG (Shanghai) Automation Equipment Co., Ltd.

Section A2,2/F, Building 6,
No. 473, West Fute 1st Road,
Waigaoqiao F.T.Z, Shanghai, 200131,
P.R. China

Phone +86 21 50375915
Fax +86 21 50375713
MP +86 15 80 07 50 53 3

xc.zeng@herzog-automation.com.cn
www.herzog-automation.com.cn

HERZOG